

Revision History Winlog.basic

Version:	Release Date
3.1.1	07.11.2024
Changed	Updated Winlog.basic to hide the internal channel, pressure unit options
Fixed	Resolved an issue where repeated use of "optimize view with Limits" caused charts to shrink by ensuring that chart range calculations account solely for measurements and limits, excluding the current y-axis range.
Changed	Removed TCP settings
Changed	Updated the system to reset the zoom level to its previous state after report generation.
Fixed	Reordered the start and stop communication threads to prevent the device search window from appearing without progressing when opening an edf2 file via command line.
Fixed	Addressed an issue where the Winlog.updater was called multiple times upon startup
Fixed	Resolved an issue where a delayed start caused a resource conflict by ensuring timers and threads start only after full initialization.
Fixed	Resolved an issue where the application sometimes crashed during measurement fixing graph moduls formatting for timespans exceeding one day.
Changed	Updated reports to preselect number format, date format, and language based on system and application settings, with options to change them before report generation.
Fixed	Corrected an issue where formatting limits in the program limit tab failed due to incorrect type access.

Version:	Release Date
3.1	23.09.2024
Fixed	Hiding one channel of a logger no longer hides the complete logger from reports
Fixed	Winlog.pro displaying Winlog.updater multiple times upon startup by increasing the default interval for update checks.
Improved	Enhanced precision of ADValue Calculation for TriggerValues and robustness for incorrectly configured loggers.
Fixed	Issue of temperature display in other precision after saving advanced event programming has been resolved.
Fixed	Issue of logger showing measurements with CRC after advanced programming options.
Fixed	Wrong date display in advanced option programming has been addressed.
Fixed	Corrected the issue where time changed by x hours in advance after saving advanced programming options.
Fixed	Detailed report now supports current view and optimal view chart selection.
Fixed	User information is now included in all reports.

Revision History Winlog.basic

Changed	Measurement remark moved from compact to detailed report for better clarity.
Fixed	Reports now support ms samplerate and associated change events.
Fixed	Issue where entering Authorization Code ends the application has been rectified.
Improved	Software now ensures logger has not been switched in interface before programming.
Fixed	Resolved the issue of the Report menu reopening a second time after printing and closing the menu.
Fixed	Corrected a translation error where “相对湿” was displayed incorrectly instead of “相对湿度” in generated Chinese reports.
Fixed	Corrected the channel index display in reports when a channel is hidden before printing by extending the edf to store the curve index and updating the channel index counting method.
Fixed	Amended the false chart legend icon in Compact Reports by extending the edf to store the curve index and reflecting hidden but non-internal channels.
Fixed	Resolved illegible comment sections in detailed reports by replacing the system font with Arial during edf2 loading to avoid display issues for older reports.
Changed	Updated reports to use language-dependent date and number formats instead of system settings based on the selected report language.
Improved	Enabled the visibility toggle of channels by clicking the "eye" icon in the channel table.
Fixed	Ensured the correct logger data is displayed in reports by collecting data from the currently selected document view and validating event start indexes for offsets.
Fixed	Ensured all data is read in packages by adjusting the process to handle missing records and aligning and concatenating data correctly without overlaps.
Fixed	Ensured consistent visibility of curves across all chart tabs and the channel list when selecting a single logger tab.
Changed	Updated the base temperature for A0 value calculations from 65°C to 70°C to align with new standards.
Fixed	Included missing report languages for Korean, Japanese, and Swedish in the software.
Fixed	Incorporated missing translations of menu items in reports, enhancing the user interface for multiple languages.
Fixed	Added a signature line to the audit trail for better record verification.
Fixed	Improved translations in Reports, ensuring all report elements are correctly displayed in the selected language.
Improved	Updated the callback level for read progress to prevent flickering of the progress bar when reading data from a full EBI 12.
Fixed	Incorrect time shifts occurring when exporting data to XLX or XLSX formats, by ensuring timestamps are correctly localized before export.
Fixed	Addressed an error occurrence when adding the Absolute Humidity Channel in Logger Channels, which also affected the Temperature Channel.

Revision History Winlog.basic

Fixed	Restored the visibility of the password reset logo in the login prompt.
Fixed	An error message appearing when saving printer changes during the report printing process has been resolved.
Fixed	Audit Trails deactivation and reactivation events not being recorded in the Reports have been addressed by adding events for activating or disabling the audit trail in Winlog.pro.
Fixed	Programming using templates not starting the logger correctly has been addressed.
Improved	Enforcing password length checks according to security settings on each login.
Improved	User calibration now takes priority over facility calibration if it is more recent, with correct localization of calibration dates.
Fixed	Display of "00:00:00" in measurement table within the Winlog.pro has been forced to ensure it is shown correctly.
Fixed	Missing Range info in detailed reports has been addressed.
Improved	Sorting by channel name now meets user expectations by forcing two-digit formatting for logger numbers.
Improved	Violation duration calculations and display in Winlog to ensure accurate measurement durations.
Fixed	The inability to change the report logo has been resolved by saving the logo setting to the registry and using a standard control for logo changes.
Fixed	Resolved issues with the Authorization Code being marked invalid during the password reset process.

Version:	Release Date
3.0	25.03.2024
Improved	Tabular report with statistics has new layout.
Changed	New reporting engine
Changed	Updated XMLParser. Now uses current Windows service
Fixed	Fileformat corruption on save has been fixed. Files from previous Winlog.pro that have this issue can be opened in the current Version.
Changed	New Setupengine and fixed setup routine
Fixed	Reinstated support for silent setup.
Improved	New device drivers for improved compatibility with Windows 10 and Windows 11.
Fixed	Inclusion of images in detailed reports has been fixed.
Improved	Userinterface now uses the latest native controls of the operating system.
Fixed	Reading issue of EBI 3x0 loggers

Revision History Winlog.basic

Fixed	Memoryleaks and crashes
Improved	Foodlogger now use two decimal places.
Fixed	Removed the display of invalid calibration dates

Version:	Release Date
----------	--------------

2.81	12.02.2019
Fixed	Repeated data in tabular reports
Fixed	Generation of tabular reports showing unwanted misleading popups

Version:	Release Date
----------	--------------

2.8	09.02.2018
Added	Support for new EBI 12 loggertypes
Improved	New PDF calibration reports
Fixed	Fixed bug when exporting data to excel

Version:	Release Date
----------	--------------

2.7	01.12.2016
Improved	Support for 1 or 2 channel TPX 310
Fixed	Fixed wrong display of chart markers when printing black/white
Changed	Removed File Explorer. Windows 10 doesn't support this control anymore
Fixed	Fixed wrong display of calibration date when using EBI 300
Fixed	Fixed wrong calibration when using EBI 300/310 with external probe activated
Added	Added Excel Export Options

Version:	Release Date
----------	--------------

2.68	06.10.2015
Improved	Improved support for core-temperature enabled loggers (EBI3x0)
Added	Added support for NFC-enabled loggers
Fixed	Fixed minor bugs

Revision History Winlog.basic

Version:	Release Date
2.67	30.07.2015
Improved	Support for EBI310 Core-Temperature-Simulation Loggers

Version:	Release Date
2.66	29.10.2014
Improved	Fixed small bug in gimp.dll introduced in 2.65

Version:	Release Date
2.65	06.10.2014
Fixed	Fixed minor bugs
Added	New measuring mode for EBI 310: Maximum measurement time
Improved	Support for external probes: TPH 500 TPH 400 TPX 220 TPX 250 TPX 310 TPC 300

Version:	Release Date
2.64	20.03.2014
Added	Added automatic software update feature
Improved	Added Support for EBI300/310 external probes

Version:	Release Date
2.63	11.07.2013
Fixed	EBI 300 and 330 loggers are programmed with 1 sec. Measurement interval when using the calibration plugin.
Fixed	English version does not display charts
Fixed	The report selection dialog shows disabled radio-buttons even if a report has been selected

Revision History Winlog.basic

Fixed	Programming EBI 330 (V1.0) with calibration plugin results in a measurement interval of 3 sec. And 100 msec.
Improved	Missing image on battery exchange wizard.
Improved	Software cuts channelnames larger than 32 characters without notification

Version:	Release Date
2.62	08.05.2013
Added	Added support for EBI20-TH-EXT
Added	Added support for MS Excel XLSX files.
Improved	Changed Chart-Behaviour: Chart axis no longer show negative values for pressure-axis and humidity axis. Furthermore, humidity axis is limited to 100%
Improved	Calibration scripts revised
Improved	Fixed wrong System-Script for setting the EBI2-Alarm-Delay
Fixed	Fixed small EBI-11 pressure-calculation issue (rounding error)
Improved	Redesigned-Reports (Tabular and Compact)
Added	Added support for sensor-serial, sensor-calibration and sensor-cycle-counter
Added	Added Support for EBI 330-T30
Added	Added Support for EBI 330-T85
Added	Added support for EBI10-TH

Version:	Release Date
2.61	14.11.2012
Fixed	User can change the tab-order of the program dialog. TC W112012264
Added	New scripts for entering Truck-Texts and alarm-delay (EBI2 Loggers)
Fixed	Different MKT-Scripts in german and english.TC W112012307
Fixed	Program dialog --> Configuration --> Mark multiple configurations Error: Button "Apply" remains active, but should be disabled in case of multiple configurations.TC W112012270
Fixed	Program dialog -> Configurations --> Select multiple configurations and press delete Error: Only the first selected configuration will be deleted TC W112012269

Revision History Winlog.basic

Fixed	- open security settings - enter invalid value for "max invalid logons" (>6) - a warning message appears indicating a value between 1 and 6 - enter a valid value and click on OK - the warning message still appears and the wrong input-field is focused TC WI12012329
Changed	Set activation energy of MKT to 83.144720. TC WI12012276
Fixed	Combobox containing measurement types is too small when using french. TC WI12012304
Fixed	Software crashes when calculating theoretical steam temperature. TC WI12012317
Improved	1st and 2nd signature can be entered independently.
Added	Added build in file explorer.
Fixed	Fixed bug when saving a file from logger that contains a user calibration
Fixed	The x-axis shows markers on the screen as well as on the report. TC WI12012311

Version:	Release Date
2.6	07.03.2012
Added	Support for EBI310 dataloggers
Added	New compact-report with additional statistics
Fixed	Limits were not used when importing Winlog 2000 files

Version:	Release Date
2.52	19.12.2011
Improved	Supported languages: German Englisch French Czech Italian Spanish Japanese Chinese Swedish
Fixed	Fixed setup bug: ReportLanguage files were missing

Version:	Release Date
2.5	18.10.2011
Added	Initial version

Revision History Winlog.basic

Version:	Release Date
1.0	20.03.2008
Fixed	Minor bugs and issues
Fixed	Issue that prevented the reset password function to work in service
Improved	More robust communication with EBI 10 / EBI 12 -Series
Improved	Compatibility with Windows 10 and Windows 11
Improved	Winlog.pro now accounts for 1000 heat cycles of EBI 12.